

Annealing Protocol

V.1 NDJ 1-22-04

1. Place the following in to microfuge tube:
 - 10ul 100uM Oligo 1
 - 10ul 100uM Oligo 2
 - 40ul H₂O
2. Boil 500ml beaker 2/3 full of H₂O.
3. Put boiling beaker in ice bucket. Put oligos in water bath.
4. Put lid on bucket and let sit for 5 minutes (melt).
5. Remove lid and let sit for 30-60 minutes (until < 40°C).
6. Dilute and use.