

Stripping Southern or Northern Blot

NDJ V.1 9.30.04

1. Prepare stripping solution: 10mM Tris pH 7.5 1%SDS

5ml	1 M Tris
5ml	10% SDS
490ml	H2O

2. Boil solution 1 liter flask (SDS will boil violently) on flame.
3. Add solution pour onto blot and cover with Styrofoam or glass plate.
4. Rock 10 minutes
5. Pour off solution
6. Repeat steps 1-5
7. Pre-hybridize and follow standard probing procedures.